


Centrifuge Instructions

- 1) Place samples to be centrifuged in the plastic inserts for the appropriate size Falcon tube (15ml on the left, 50 ml on the right).

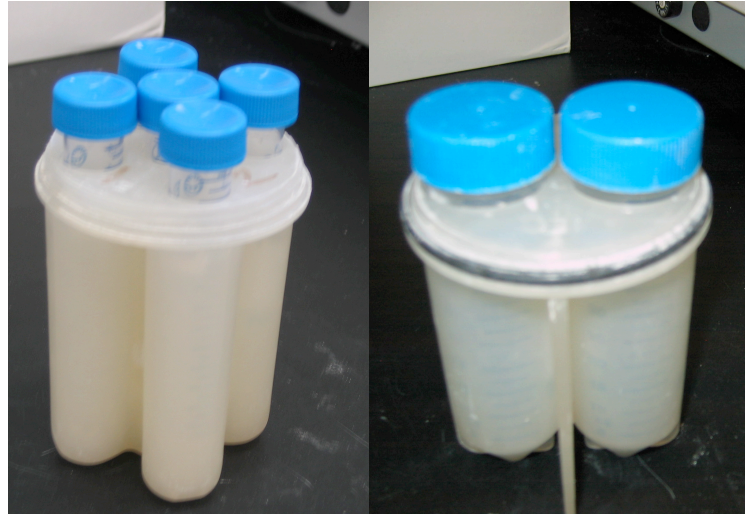


- 2) Turn on the centrifuge.
- 3) Press the release button to open the lid. The power must be on to open the lid.



- 4) Place the inserts into the centrifuge cups. **MAKE SURE THAT THE CENTRIFUE IS EQUALLY WEIGHTED ON ALL SIDES.** If the centrifuge is not balanced, it will shut itself down.



- 5) Set the centrifuge to the desired rotations per minute (RPM) and length of time to be centrifuged. For carbonate sample 1 minute at 4000 RPMs is generally adequate. **Do not adjust the rotor/radius.**

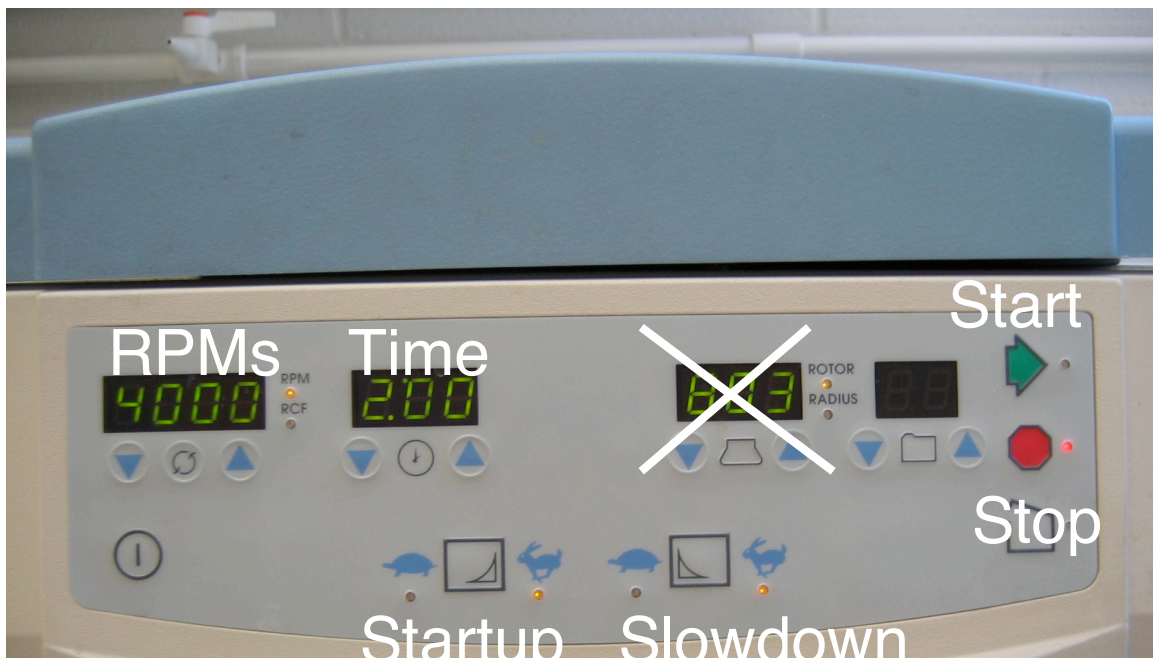

- 6) Make sure that both the startup and slowdown settings are set to fast (the rabbit icon).
- 7) Press the green arrow to start
- 8) If you need to stop the run before the time is finished, press the red stop sign button to abort.
- 9) Press the release button to release the lid.